

Summary: Prior to analysis, samples are digested via Kjeldahl digestion to hydrolyze phosphorus to orthophosphate. Orthophosphate reacts with molybdenum(VI) and antimony(III) in an acidic medium to form an antimony-phosphomolybdate complex. This complex is subsequently reduced with ascorbic acid to form a blue color, and the absorbance is measured at 880 nm.

Interferences: The presence of more than 50 mg/L ferric iron, more than 10 mg/L copper, or more than 10 mg/L silica may interfere. Samples with background absorbance at the analytical wavelength may interfere. Filter turbid samples prior to analysis.


Performance Specifications:

Range:	0.02–20 mg/L P
Throughput:	48 samples/hour
Precision:	
0.10 mg/L	<1% RSD
4.0 mg/L	<1% RSD
Method Detection Limit (MDL):	0.005 mg/L P

Chemicals:

Ammonium Molybdate Tetrahydrate, (NH ₄) ₆ Mo ₇ O ₂₄ •4H ₂ O	Hydrochloric Acid, concentrated, HCl
Antimony Potassium Tartrate Hemihydrate, K(SbO)C ₄ H ₄ O ₆ •½H ₂ O	Potassium Phosphate Monobasic, KH ₂ PO ₄
Ascorbic Acid, C ₆ H ₈ O ₆	Potassium Sulfate, K ₂ SO ₄
Chloroform, CHCl ₃	Red Mercuric Oxide, HgO
DOWFAX® 2A1 (OI Analytical Part #A000080)	Sodium Chloride, NaCl
	Sodium Hydroxide, NaOH
	Sulfuric Acid, concentrated, H ₂ SO ₄

Basic Flow Diagram:

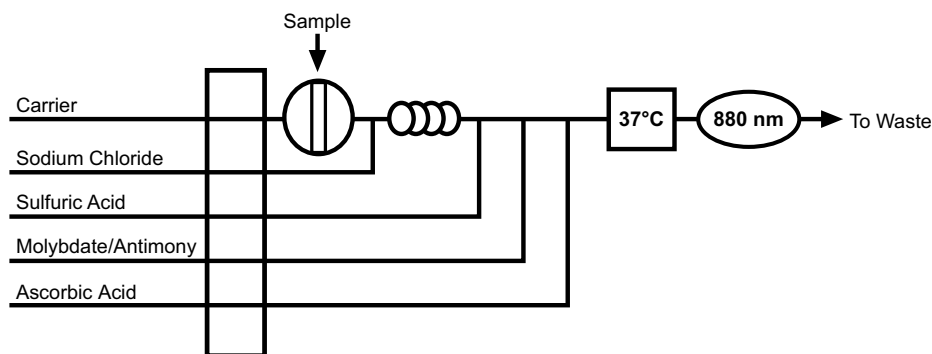

Note: This method complies with USEPA Method 365.1

Selected References: *Methods for Chemical Analysis of Water and Wastewater*; EPA-600/4-79-020; U.S. Environmental Protection Agency, Office of Research and Development, Environmental Monitoring and Support Laboratory: Cincinnati, OH, 1984; Method 365.1.

Standard Methods for the Examination of Water and Wastewater, 20th ed.; American Public Health Association: Washington, D.C., 1998.

DOWFAX is a registered trademark of Dow Chemical.