

Summary: Prior to analysis, samples are digested via Kjeldahl digestion to hydrolyze phosphorus to orthophosphate. Orthophosphate reacts with molybdenum(VI) and antimony(III) in an acidic solution to form an antimony-phosphomolybdate complex. This complex is subsequently reduced with ascorbic acid to form a blue color, and the absorbance is measured at 880 nm.

Interferences: Filter turbid samples prior to analysis. Samples with background absorbance at the analytical wavelength may interfere.


Performance Specifications:

Range:	0.01–5.0 mg/L
Throughput:	45 samples/hour
Precision:	
0.10 mg/L	<5% RSD
1.0 mg/L	<1% RSD
Method Detection Limit (MDL):	0.002 mg/L

Chemicals:

Ammonium Molybdate Tetrahydrate, (NH ₄) ₆ Mo ₇ O ₂₄ •4H ₂ O	Hydrochloric Acid, concentrated, HCl
Antimony Potassium Tartrate Hemihydrate, K(SbO)C ₄ H ₄ O ₆ •½H ₂ O	Potassium Phosphate Monobasic, KH ₂ PO ₄
Ascorbic Acid, C ₆ H ₈ O ₆	Potassium Sulfate, K ₂ SO ₄
DOWFAX® 2A1 (OI Analytical Part #A000080)	Red Mercuric Oxide, HgO
	Sodium Hydroxide, NaOH
	Sulfuric Acid, concentrated, H ₂ SO ₄

Basic Flow Diagram:

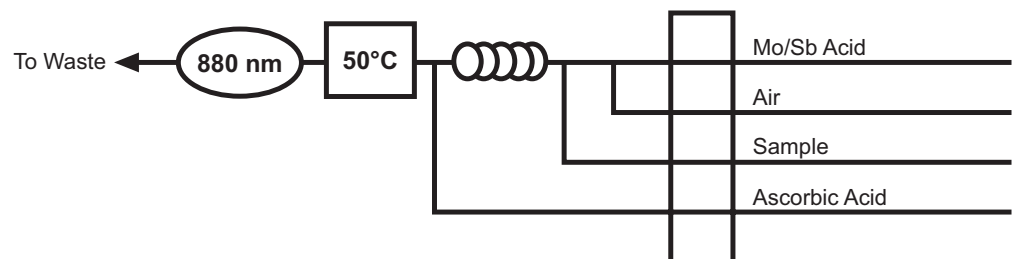

Note: This method complies with USEPA Method 365.1.

Selected References: *Methods for Chemical Analysis of Water and Wastewater*; EPA-600/4-79-020; U.S. Environmental Protection Agency, Office of Research and Development, Environmental Monitoring and Support Laboratory: Cincinnati, OH, 1984; Method 365.1.

Standard Methods for the Examination of Water and Wastewater, 20th ed.; American Public Health Association: Washington, D.C., 1998.

DOWFAX is a registered trademark of Dow Chemical.